

***Toughest Samples.
Fastest Results.***

OMNI BEAD RUPTOR

Bead Mill Homogenizer



- Convenient Front-Loading Design
- Consistent & High-Quality Results
- No Cool-Down Between Runs
- Broadest Performance Range: 0.8m/s—8m/s
- Compatible with 0.5mL, 2mL & 7mL Sample Tubes



www.omni-inc.com

The **Omni Bead Ruptor 24** is the newest generation of bead mill homogenizers. This system is specifically designed for the grinding, lysing and homogenization of biological samples prior to any molecular extractions.

Using sample tubes pre-filled with a variety of lysing beads, the Bead Ruptor 24 vigorously and uniformly shakes the tubes providing an efficient, consistent, high-yield and quality homogenization usually in less than 40 seconds.

The Omni Bead Ruptor 24 is an ideal solution for releasing DNA, RNA, proteins, enzymes, etc. from very tough samples while still retaining molecular integrity.



Product Features:

- Simultaneously homogenize either 24 x 2mL or 12 x 7mL tubes
- Convenient front-loading tube holder position for ease of use and optimal bead/sample interactions
- Broad speed range for any type of sample processing
- High reproducibility: number of cycles, time, speed, dwell/pause
- No Cool-down required between runs: process hundreds of samples per day
- 99 programmable memory settings
- Advanced sample temperature monitoring system
- Multi-lingual user interface
- Disposable tubes ensure no threat of cross-contamination or sample degradation

Omni Bead Ruptor Lysing Tubes are available pre-filled with a variety of bead types.

Our lysing tubes are ideally suited for nucleic acid, drug, proteins, enzymes, virus etc. extractions. **Tubes are RNase & DNase free.**

Efficiently lyse with or without buffer:

- Human
- Animal
- Plant and food
- Micro-organisms
- Soil and sediments
- Forensic, swabs, feces
- Drugs, polymers, lotions etc.



1.4mm & 2.8mm Ceramic Beads

Ideal for soft animal, plant or human tissues:

- brain
- liver
- skin
- leaves
- gonad
- muscle, etc.



0.15mm & 0.7mm Garnet Beads

"Universal" bead mix

- Good for tough samples
- Not for use with soil



2.38mm Metal Beads

Ideal for any hard tissue types:

- hair
- teeth
- bones
- seeds
- corn
- rice



0.25 Carbide Beads

Ideal for microbial and soil RNA



0.1mm & 0.5mm Glass Beads

Ideal for microorganisms

Optimal for:

- spores
- fungi
- yeast
- e-coli, etc.




Unique Bead Ruptor Tube Rack for sample storage down to -80°C.

The Omni Bead Ruptor 24 is ideally suited for:

- DNA / RNA electrophoresis
- Northern, Southern, Western blot analysis
- Liquid chromatography
- Enzyme isolation
- Protein purification and production
- Food safety, GMO detection
- Cosmetic preparation
- Environmental surveys
- BSE Screenings
- Drug detections
- qPCR, RT-PCR
- And many more....



Before & After

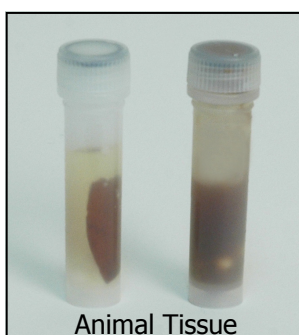

Specifications:

Processing Range: 25 μ L to 5mL

Power Rating: 100-230 VAC/ 50-60 Hz

Dimensions: 15.5" (W) x 17" (D) x 13" (H) [24" tall with lid open]

Weight: 24 kg (54 lbs.)

Compliance: CE Compliance

Run Time: 1 second—9:59 minutes

Number of Programs: 99 Saved Programs

Number of Cycles: 1—10

Performance Range: 0.8m/s—8m/s by 0.15m/s increments

Below are recommended settings for the Bead Ruptor 24.

Actual settings will vary based on sample conditions.

Detailed application notes can be found at

www.omni-inc.com/applications

		BEAD MATRIX	SETTING	APPLICATION
Sample Name and Type	Lysing matrix	Speed – number of cycle x time - pause	Note	
MICROBIAL	Bacteria	Garnet 0.15mm	3.25m/s – 1 x 5 min – 0s	DNA isolation
ANIMAL	Muscle	Ceramic 2.8mm	5.65m/s – 2 x 45 sec. – 30 sec.	RNA isolation
	Brain	Ceramic 2.8mm	5.65m/s – 2 x 45 sec. – 30 sec.	RNA isolation
	Liver	Ceramic 2.8mm	5.65m/s – 2 x 45 sec. – 30 sec.	RNA isolation
	Kidney	Ceramic 2.8mm	5.65m/s – 2 x 45 sec. – 30 sec.	RNA isolation
	Lung	Ceramic 2.8mm	5.65m/s – 2 x 45 sec. – 30 sec.	RNA isolation
	Heart	Ceramic 2.8mm	5.65m/s – 2 x 45 sec. – 30 sec.	RNA isolation
	Spleen	Ceramic 2.8mm	5.65m/s – 2 x 45 sec. – 30 sec.	RNA isolation
	Leaves	Ceramic 2.8mm	6.95m/s – 2 x 45 sec. – 30 sec.	RNA isolation
PLANT	Leaves	Metal 2.38mm	3.25m/s – 1 x 3 min	DNA isolation
	Seeds	Metal 2.38mm	4.7m/s – 1 x 3 min	DNA isolation
	Stems	Ceramic 2.8mm	6.95m/s – 2 x 45 sec. – 30 sec.	RNA isolation
	Stems	Metal 2.38mm	3.55m/s – 1 x 3 min.	DNA Isolation
	Roots	Ceramic 2.8mm	6.95m/s – 2 x 45 sec. – 30 sec.	RNA isolation
	Root	Metal 2.38mm	4m/s – 1 x 3 min	DNA isolation
	Flowers	Ceramic 2.8mm	6.95m/s – 2 x 45 sec. – 30 sec.	RNA isolation
	Soil	Garnet 0.7mm	3.25m/s – 1 x 5 min	DNA isolation



935-C Cobb Place Blvd.
 Kennesaw, GA 30144
 Phone: 800-776-4431
 Fax: 770-421-0206
omni@omni-inc.com

www.omni-inc.com